

Document Title: Standards Insert, Hang Tag		Part # and Rev. 18029	
		Release Date: 11/06/15	
Rev.	Description	Revision By	Date
Date -- Time	Uploaded Quick Start to repository at http://publications.decagon.com . Please ask archivist for previous versions.	NB	11/06/15

Production Filename:

http://publications.decagon.com/Inserts/18029_Standards_Insert_Hang_Tag-4up_Print.pdf

Dimensions: 2.25 inch wide, 6 inch tall

Colors: Full Color

Materials: 10223

Finish: N/A

Bindery: Drill holes, Trim, Finish

Packaging: N/A

Special Notes: 1” Hole 0.25” from edge with the center of the hole at y:0.75” x:1.125”. Fold at 1.75” from top edge.

6Z081	51-11-90	<p>Instructions Read First</p> <ol style="list-style-type: none"> 1. Replace the flat cap with the nozzle cap and ensure it is well secured. 2. Hold the red cap and shake well. 3. Take a clean kimwipe and wipe up the nozzle as you remove the red cap from the nozzle tip. 4. Drip two or three drops into a waste container. Make sure not to touch the nozzle tip to the container. 5. Fill a sample cup until the bottom of the cup is covered. 6. Recap the standard bottle. Put the lid on the sample cup and label it. Under most conditions, the sample cup containing the standard can be used for up to 4 hours if kept covered and sealed with parafilm when not being read. 7. Make sure the red cap remains on the standard bottle at all times when not in use. 8. There might be some salt crystallization around the cap and nozzle for .760_{aw} standard. It can be wiped off with a tissue and disposed of in the trash. Do not rinse the bottle. 	<p>Have Questions? ☎ 1-509-332-5601 ✉ support@aqualab.com</p>
-------	----------	---	---